

MetAmino Kit®

Installation manual for GC-MS MetAmino® kit



Installation Manual

Ver. EN221111



©Copyright Chromservis s.r.o. All rights reserved.

Important notes: **Read this manual carefully before the installation.**

1. Contents of the Manual

1. Contents of the Manual.....2

2. Introduction3

3. Installation.....3

4. Manufacturer Contact4


2. Introduction

The ZIP file includes MS spectra of MetAmino[®] derivatives used in the GC-MS kit. You can download this file at <https://www.chromservis.eu/en/metamino-gc-ms-nist-library>.


3. Installation

- Please unzip the file MetAminoGC.
- Copy the unzip file to the NISTMS folder, typically: ProgramFiles/NISTMS/MSSEARCH".Extended box (reagents, liquid media, chemicals and accessories for an additional 300 samples)


•

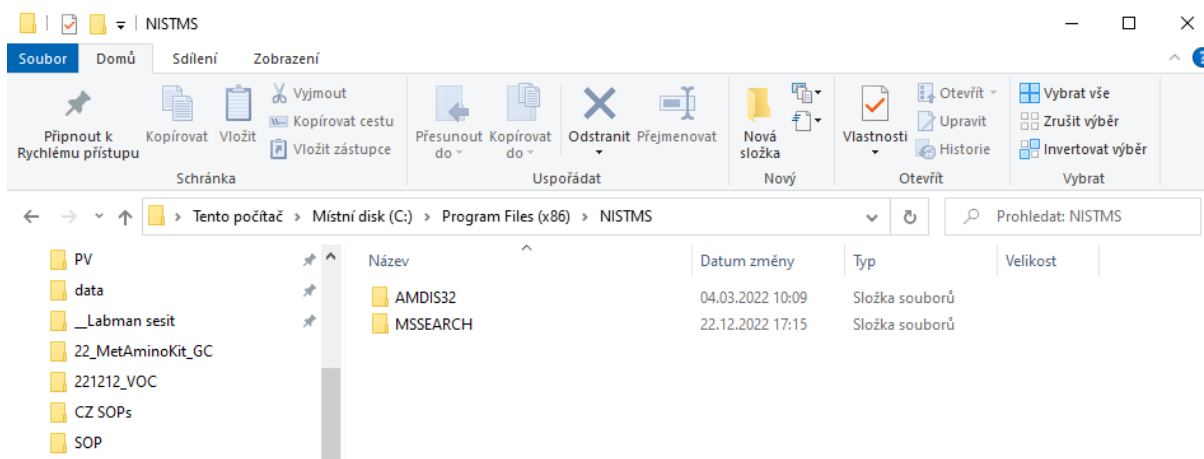

- Open the program "NIST MS Search".
- In main menu of NIST MS Search click on the Options and then click on Library Search Options.

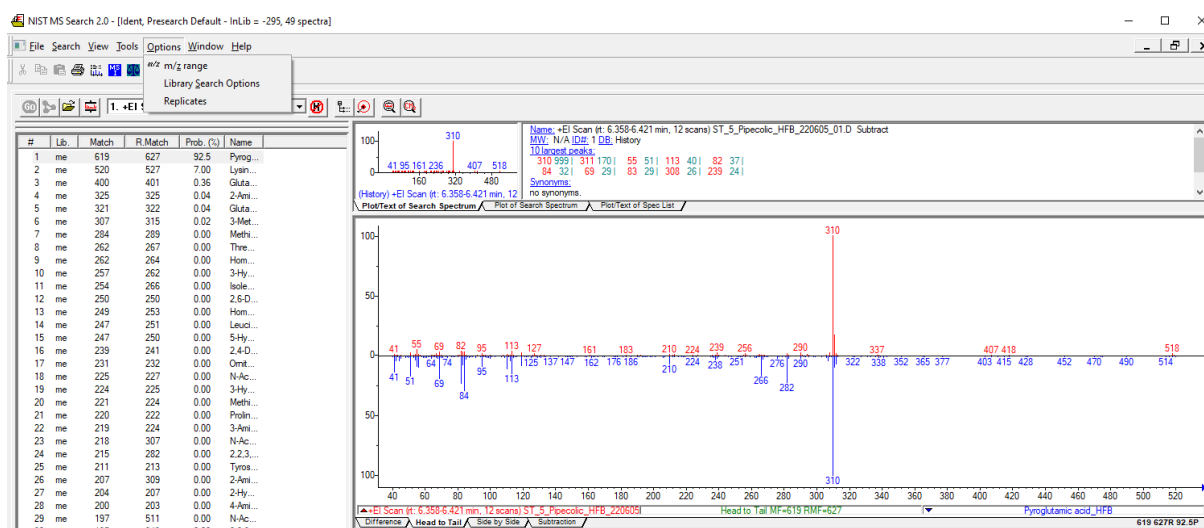

- In the window Library Search Options select the „metaminogc” and press the button Add for include the metaminogc library.

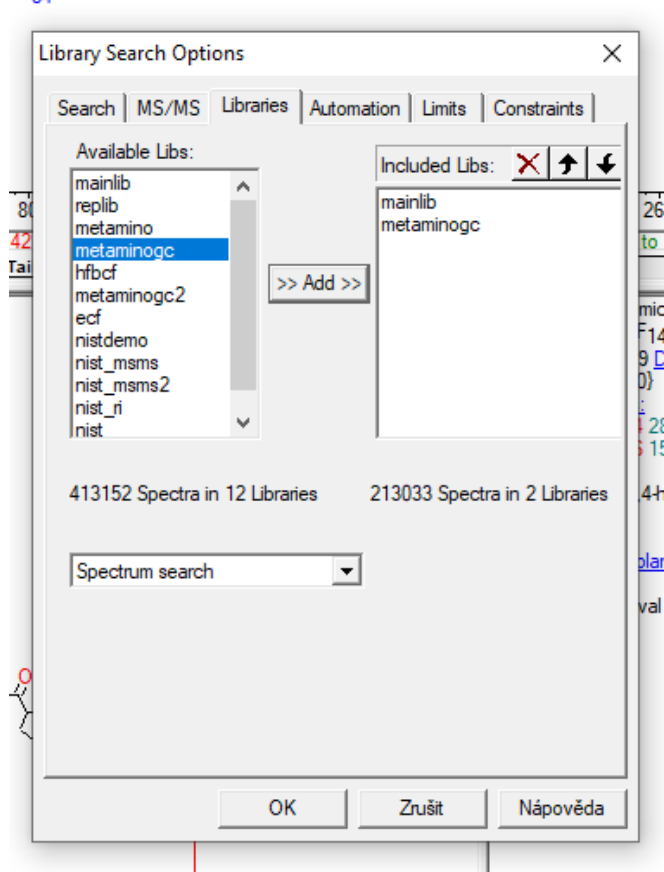

4. Manufacturer Contact

Chromsevis s.r.o.
 Jakobiho 327
 109 00 Praha 10 – Petrovice
 The Czech Republic
www.chromservis.eu

The MetAmino® is a trademark registered by Chromservis s.r.o. Other trademarks and trade names may be used in this document to refer to either the entities claiming the marks and/or names or their products and are the property of their respective owners. We disclaim proprietary interest in the marks and names of others.

No part of this document may be copied, translated or changes without the written consent of Chromservis s.r.o. Chromservis s.r.o. reserves the rights of change or improvement of this User manual.